
Precision Fresnel Biprism Interference Apparatus

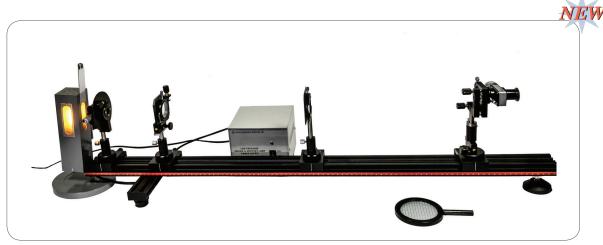



- Perform Fresnel's Biprism Experiment on a 1m optical bench
- Precision optics allow fringe position measurements to 1/100mm
- Easy alignment and adjustment of the two virtual light sources

Fresnel's Biprism Experiment is the fundamental example of common path interference by division of wavefront. Light from a slit passes through side-by-side prisms with very shallow angles to separate it into two almost parallel beams. When these beams are brought to overlapping by a convex lens, they interact as if they originated from two virtual sources side by side and generate equally-spaced fringes. The original version of the experiment by Fresnel is difficult to perform because it requires precise adjustment of the slit and prisms, a dark room, and a large amount of space to separate the fringes for easy viewing.

This apparatus allows these issues to be resolved by using a precision angle-adjustable slit, a prism carrier with both lateral and angular adjustment, and a micrometer eyepiece to allow fringe separation measurements to 1/100mm. The experiment can be carried out on a 1m optical bench with only dimmed lighting.

The micrometer eyepiece has a range of ±4.00 mm with a backlash-free, easy-to-read drum scale.

During setup, the two virtual sources are focused onto a ground glass screen for easy alignment and adjustment so that the sources are parallel and of equal intensity.

The fringes show good contrast and separation for dark-dark position measurements.

Contents:

1. Optical Bench, 1000mm	ı
2. Riders, Lateral Adjustment	2
3. Riders, Plain	2
4. Sodium Lamp, 20W, 110V	
5. Adjustable Slit	
6. Biprism Carrier, Adjustable	
7. Lens Carrier, Plain	
8. Eyepiece Carrier, Adjustable	
9. Micrometer Eyepiece	
10. Ground Glass Screen	
11. Eyepiece Adapter Ring	
12. Height Fixing Collar 3	3
13. Biprism	
14. Lens. f = +150mm	ı

 Item No.
 Description

 PFB001
 Precision Fresnel Biprism Interference Apparatus